




DIAPHRAGM VACUUM PUMPS

DATA SHEET E 040



N 860 FTE

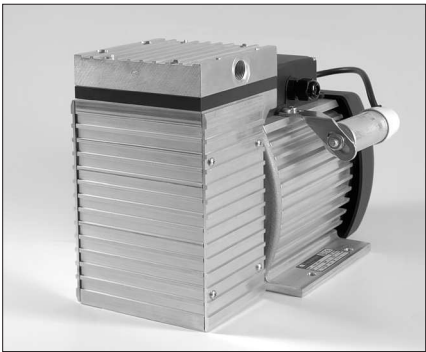

N 860 ANE



N 860.3 ANE

Concept

The Diaphragm Vacuum Pumps from KNF are based on a simple principal - an elastic diaphragm, fixed on its edge, moves up and down its central point by means of an eccentric. in this way the substance is transferred using automatic valves.

The pumps are equipped with the patented stress-optimised structured diaphragm, resulting in a high pneumatic performance, long product life and compact size. Special valves ensure that the product can cope easily with vapour and condensation.

Thanks to the KNF modular system, the parts used to transfer the gases can be made from materials with varying degrees of resistance. The customer has a choice of pump drives ranging from a selection of A/C motors to explosion-proof models. Please contact us for further details.

Features

- Uncontaminated flow**
No contamination of the media due to oil-free operation
- Maintenance-free**
- Compact size**
due to structured diaphragm
- High performance**
because of structured diaphragm
- High level of gas tightness**
- Chemically resistant versions**
- Long product life**
thanks to structured diaphragm
- Very quiet and little vibration**
- Copes well with vapour and condensation**
- Starts against vacuum**
- Cool running motor**
even when in constant use
- Can operate in any installed position**

Areas of use

The Diaphragm Vacuum Pumps offer a high level of performance despite their small size, as well as an excellent price performance ratio. They are required especially in the fields of analysis, medicine and production technology.

The pumps are used for sucking gases, taking samples (even liquids in a vacuum) and evacuating vessels.

PERFORMANCE DATA

Type	Delivery (l/min)	Vacuum (mbar abs.)	atm. Press.	Pressure (bar g)	Weight (kg)
N 860 ANE	60	80		1	12.7
N 860 ATE	60	80		1	12.7
N 860 FTE	60	80		1	12.5
N 860.3 ANE	60	2		1	14.7
N 860.3 ATE	60	2		1	14.7
N 860.3 FTE	60	2		1	14.3

N 860 ANE N 860 ATE

PERFORMANCE DATA

Type and Order No.	Delivery at atm. Pressure (l/min) ¹⁾	Max. operating pressure (bar g)	Ultimate vacuum (mbar abs.)
N 860 ANE	60	1	80
N 860 ATE	60	1	80

¹⁾ Litre at STP


MOTOR DATA

Protection class	IP 54		
Voltage/Frequencies (V/Hz)	~230/50		
Power P ₁ (W)	200		
Operating current (A)	1.5		


MODEL CODES AND MATERIALS

Type and Order No. ⁴⁾	Pump head	Diaphragm	Valves
N 860 ANE	Aluminium	EPDM	EPDM
For slightly aggressive or corrosive gases and vapours			
N 860 ATE	Aluminium	PTFE coated	FFPM

²⁾ See also „MODEL CODE FOR EASY ORDERING“

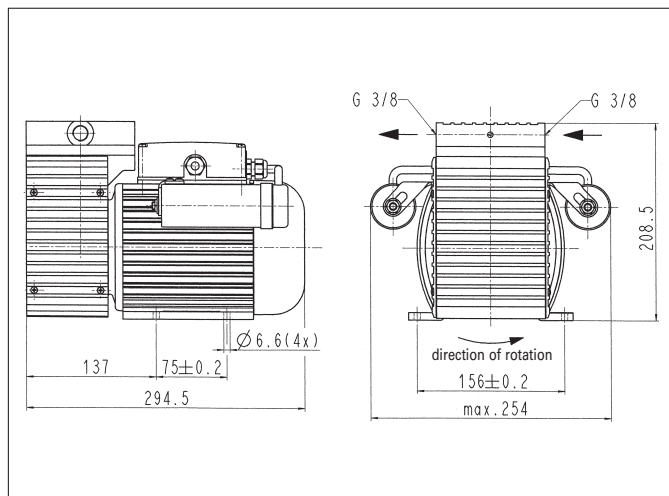Dimensions ³⁾ (mm)

³⁾ All dimensional tolerances conform to DIN ISO 2768-1, Tolerance Class V

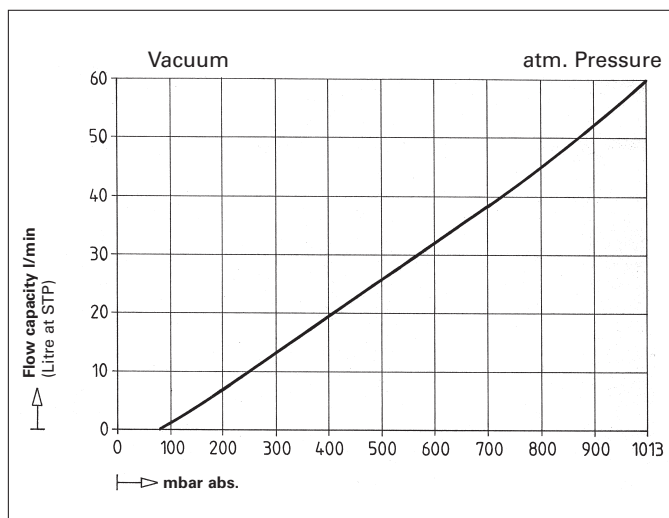

N 860 FTE

Chemically resistant version

PERFORMANCE DATA

Type and Order No. **	Delivery at atm. pressure (l/min) ¹⁾	Max. operating pressure (bar g)	Ultimate vacuum (mbar abs.)
N 860 FTE	60	1	80



¹⁾ Litre at STP

MOTOR DATA

Protection class	IP 54		
Voltage/Frequencies (V/Hz)	~230/50		
PowerP ₁ (W)	200		
Operating current (A)	1.5		

MODEL CODES AND MATERIALS

Type and Order No. ²⁾	Pump head	Diaphragm	Valves
N 860 FTE	PTFE	PTFE coated	FFPM

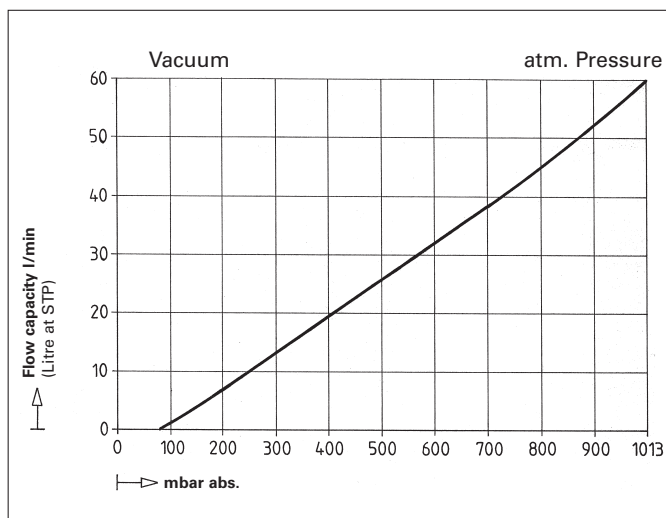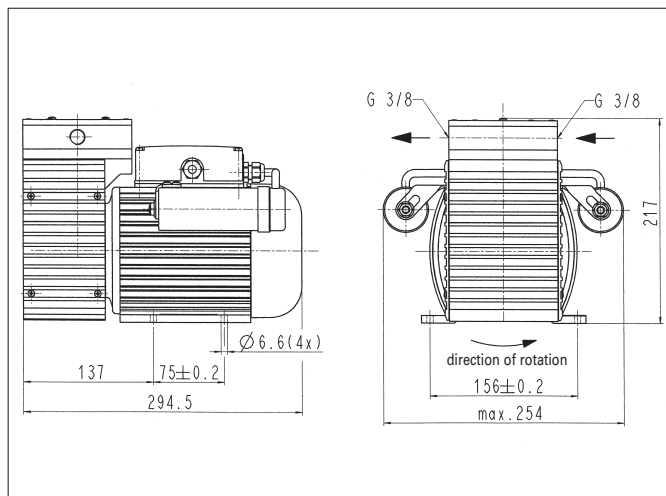

N 860.3 ANE N 860.3 ATE

N 860.3 FTE

Chemically resistant version

Type and Order No.	Delivery at atm. pressure (l/min) ¹⁾	Max. operating pressure (bar g)	Ultimate vacuum (mbar abs.)
N 860.3 ANE	60	1	2
N 860.3 ATE	60	1	2

¹⁾ Litre at STP

Protection class	IP 54		
Voltage/Frequencies (V/Hz)	~230/50		
Power P ₁ (W)	220		
Operating current (A)	1.6		





Type and Order No. ²⁾	Pump head	Diaphragm	Valves
N 860.3 ANE	Aluminium	EPDM	EPDM
For slightly aggressive or corrosive gases and vapours			
N 860.3 ATE	Aluminium	PTFE coated	FFPM

Type and Order No.	Delivery at atm. pressure (l/min) ¹⁾	Max. operating pressure (bar g)	Ultimate vacuum (mbar abs.)
N 860.3 FTE	60	1	2


¹⁾ Litre at STP

Protection class	IP 54		
Voltage/Frequencies (V/Hz)	~230/50		
Power P ₁ (W)	220		
Operating current (A)	1.6		

Type and Order No. ²⁾	Pump head	Diaphragm	Valves
N 860.3 FTE	PTFE	PTFE coated	FFPM

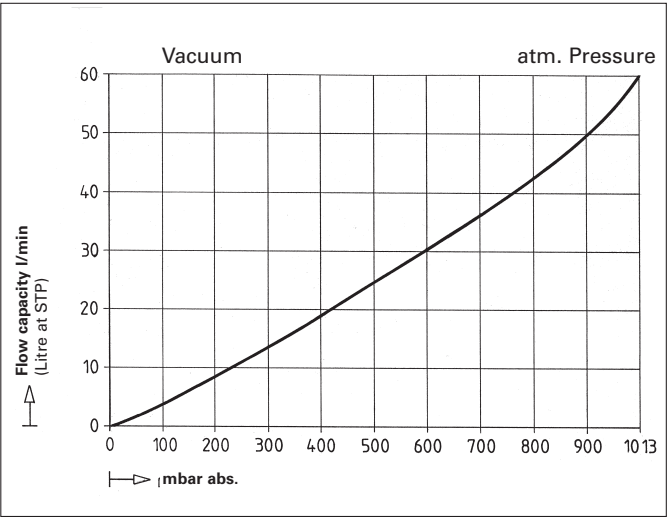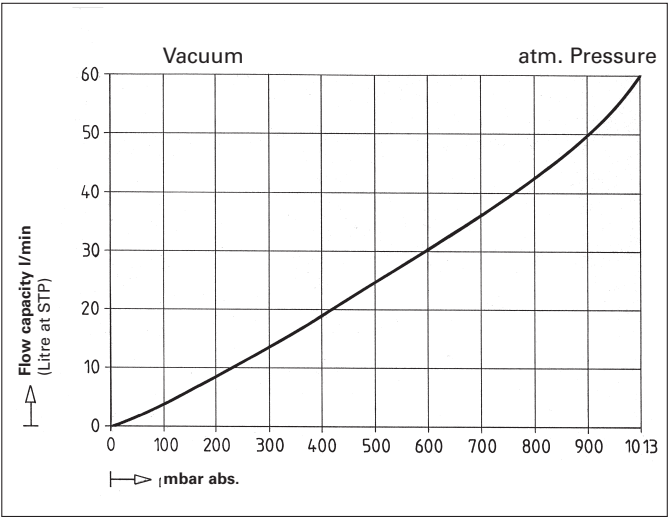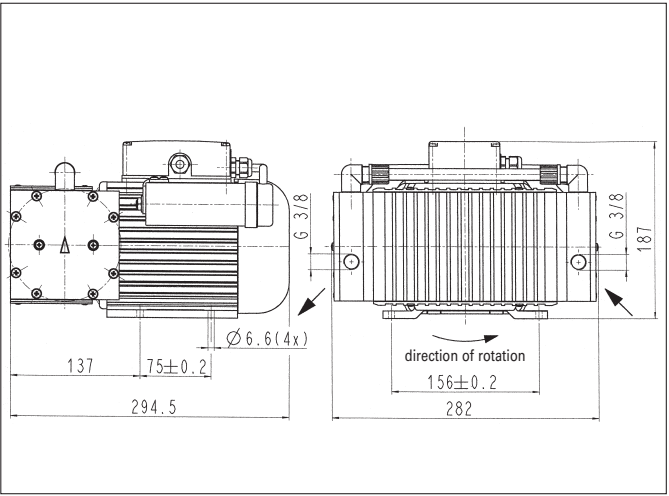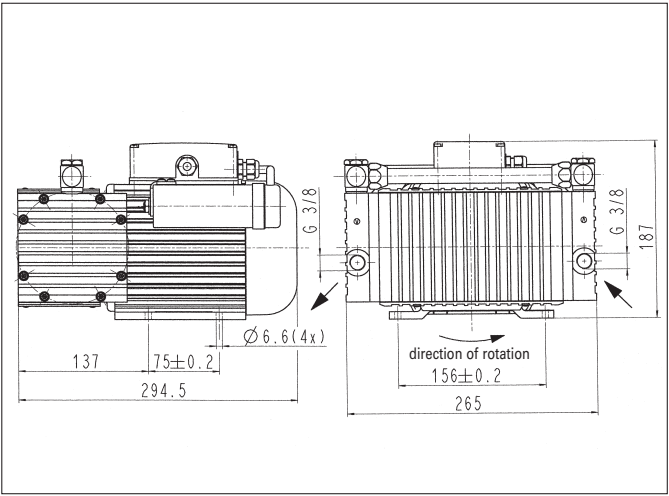


Pump down time for 20 l receiver

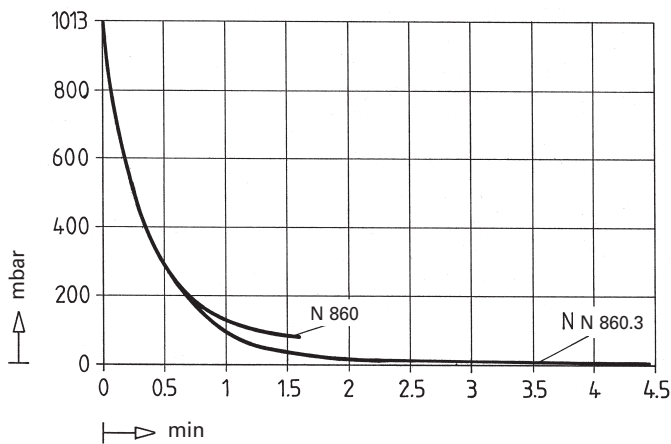

MODEL CODE FOR EASY ORDERING

The model code is identical to the order number. It is set up as follows:

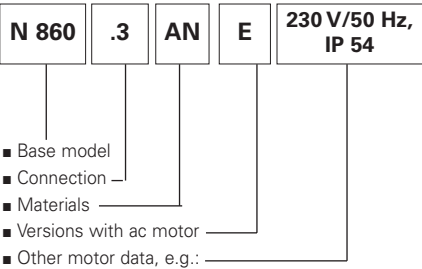

In addition the motor data must be given in the purchase order (voltage, frequency, and protection class). In our extensive program you can be sure to find the pump you need for your particular application.

Accessories

Description	Order No.	Details
Silencer	045993	G 3/8
Filter	046164	G 3/8
Hose connector	045859	G 3/8 / Ms
Hose connector	046851	G 3/8 /PVDF
Gasket	044982	

TECHNICAL DETAILS

Maximum permissible gas and ambient temperature: between +5°C and +40°C.

Motors with other voltages, frequencies and protection classes on request.

Hints on function, installation, and service: see back side.

KNF - the competent partner for vacuum and compressor technology. Especially for unusual problems. Call us and talk to our application engineers.


KNF Neuberger GmbH
Diaphragm Pumps + Systems
Alter Weg 3
D 79112 Freiburg
Tel. ++49 (0)7664/5909-0
Fax ++49 (0)7664/5909-99
E-mail: info@knf.de
www.knf.de

HINTS ON FUNCTION, INSTALLATION, AND SERVICE

FUNCTION OF KNF DIAPHRAGM VACUUM PUMPS

An elastic diaphragm is moved up and down by an eccentric (see illustration). On the down-stroke it draws the air or gas being handled through the inlet valve. On the up-stroke the diaphragm forces the medium through the exhaust valve and out of the head. The compression chamber is hermetically separated from the drive mechanism by the diaphragm. The pumps transfer, evacuate and compress completely oil-free.

Diaphragm pump

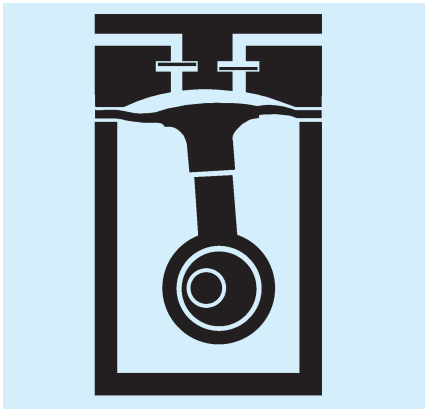

HINTS ON INSTALLATION AND OPERATION

- Range of use: Transferring air and gases at temperatures between +5°C and +40°C
- Permissible ambient temperature: between +5°C and +40°C
- The standard pumps are not suitable for use in areas where there is a risk of explosion. In these cases there are other products in the KNF program - please ask us for details
- The pumps are designed to start against vacuum. Pumps that start against pressure are available on request.
- To prevent the maximum operating pressure being exceeded, restriction or regulation of the air flow should only be carried out in the suction line
- Components connected to the pump must be designed to withstand the pneumatic performance of the pump
- Install the pump so that the fan can draw in sufficient cooling air
- Fit the pump at the highest point in the system, so that condensate cannot collect in the head of the pump - that prolongs working-life.

HINTS ON SERVICE

The diaphragm and valve plates are the only parts of the KNF diaphragm pumps subject to wear. They are easy to change, as no special tools are needed.

If you have any questions, please call our application engineers (see below for contact telephone number).

YOUR LOCAL PARTNER AROUND THE WORLD

KNF SALES ORGANIZATION

Australia KNF Regional Office

Moreland West VIC 3055
Tel. +61 3 9386 4959
info@knf.com.au
www.knf.com.au

Benelux Netherlands KNF Verder B.V.

3451 GG Vleuten
Tel. +31 30 677 92 40
info@knf-verder.nl
www.knf-verder.nl

Benelux Belgium and Luxemburg

KNF Verder N.V.
2630 Aartselaar
Tel. +32 3 871 96 24
info@knf.be
www.knf.be

China KNF Technology (Shanghai) Co., Ltd.

Shanghai 201203
Tel. +86 21 5109 9695
info@knf.com.cn
www.knf.com.cn

France KNF Neuberger SAS

68128 Village-Neuf
Tel. +33 389 70 35 00
info@knf.fr
www.knf.fr

Germany KNF Neuberger GmbH

79112 Freiburg
Tel. +49 7664 5909 0
info@knf.de
www.knf.de

India KNF Pumps + Systems (India) Pvt. Ltd.

Hinjewadi Pune 411 057
Tel. +91 20 640 13 923
info@knfpumps.in
www.knfpumps.in

Italy KNF Italia S.r.l.

20063 Cernusco s. Naviglio (MI)
Tel. +39 02 929 04 91
info@knf.it
www.knf.it

Japan KNF Japan Co. Ltd.

Tokyo 104-0033
Tel. +81 3 3551 7931
info@knf.co.jp
www.knf.co.jp

Korea KNF Neuberger Ltd.

135-502 Seoul
Tel. +82 2 959 0255
knf@knfkorea.com
www.knfkorea.com

Latin America KNF Regional Office

Tel. +1 609 649 1010
gb@knf.com
www.knf.com/es

Morocco, Tunisia, Algeria

KNF Neuberger SAS
68128 Village-Neuf
Tel. +33 389 70 35 00
info@knf.fr
www.knf.fr

Singapore KNF Regional Office

Tel. +65 9722 1994
info@knf.com.sg
www.knf.com.sg

Sweden, Finland, Denmark, Norway

KNF Neuberger AB
117 43 Stockholm
Tel. +46 8 744 51 13
info@knf.se
www.knf.se

Switzerland KNF Neuberger AG

8362 Balzerswil
Tel. +41 71 973 99 30
knf@knf.ch
www.knf.ch

UK, Ireland KNF Neuberger U.K., Ltd.

Witney, Oxfordshire OX28 4FA
Tel. +44 1993 77 83 73
info@knf.co.uk
www.knf.co.uk

USA, Canada KNF Neuberger, Inc.

Trenton, NJ 08691-1810
Tel. +1 609 890 8600
knfusa@knf.com
www.knfusa.com

KNF PRODUCT CENTERS

Gas Pumps KNF Neuberger GmbH

DE-79112 Freiburg
info@knf.de
www.knf.de

Micro Gas Pumps KNF Micro AG

CH-6260 Reiden
info@knf-micro.ch
www.knf-micro.ch

Liquid Pumps KNF Flodos AG

CH-6210 Sursee
info@knf-flodos.ch
www.knf-flodos.ch